

**DNA WI-37 Analysis of Vendor Generated Profiles per
DNA SOP 21**

Approved by Director: Dr. Guy Vallaro

Document ID: 8131
Revision: 3
Effective Date: 5/24/2019
Status: Published
Page 1 of 1

1. See SOP 21.7 for deconvolution and statistical analysis of F6C profiles. See SOP 31.10 for general protocols regarding the qualitative interpretation of vendor generated F6C profiles.
2. If necessary, the assigned DSS analyst may import the relevant sample files into GeneMarker HID. The analyst may review the analyzed data from the vendor but will not make any changes to the data.
3. The vendor lab analytical thresholds are described below for reference:

A) Bode Cellmark

Color	RFU
B	100
G	100
Y	100
R	100
P	100

B) DNA Labs International

Color	RFU
B	84
G	93
Y	39
R	42
P	64

4. Determine the number of contributors as described in SOP 31.7. Non-artifact, sub-analytical peaks are used to assess the number of contributors $\geq$ the following RFU levels for each vendor lab where the level of machine noise and the quality of the profile allows:

a) Bode Cellmark

Color	RFU
B	50
G	65
Y	50
R	50
P	50

B) DNA Labs International

Color	RFU
B	25
G	25
Y	25
R	25
P	25

5. Document Batch Paperwork review on QR-4a and Case Review on QR-4.