

Approved by Director: Dr. Guy Vallaro

A. Purpose: To establish maintenance protocols for the latent print section's Cyanosafe superglue processing chambers.

B. Responsibility:

C. Procedure:

1. For superglue chamber operation and maintenance refer to the operator's manual and maintenance log kept in the latent print processing laboratory.
2. Existing filters have an additional 30 days remaining after reaching "0" on the chamber's display screen. New filters may be ordered through the Sirchie company.
3. Replace the carbon filter as described in the Cyanosafe operator's manual. The manual is kept in the Cyanosafe processing room by the superglue control log.
4. Contact AirClean Systems (1-800-849-0472) and give control number to service attendant. (The control number will be located on the edge of the filter). The service attendant will provide a serial number, which needs to be inputted into each unit.
5. See "Directions For Entering New Serial Number" document in Operator's Manual.
6. Log information onto maintenance sheet. Chambers not operating properly will be removed from service until repaired or replaced.

D. Sources of Error:

E. References:

Cyanosafe Instruction Manual